

## Enclosed Plate and Frame Filter BECO INTEGRA® PLATE 600 EC

### Stainless Steel

BECO INTEGRA PLATE 600 EC enclosed plate and frame filter is the depth filtration system for critical filtration tasks in the pharmaceutical and biotech, chemical, cosmetics, and food industries.

BECO INTEGRA PLATE 600 EC is a multi-sheet-filter with an external chamber. The product channels are sealed with depth filter sheets only. The sealing of the whole filter package is realized by means of gaskets around the external chamber.

#### Special Features

- **CIP** of filter package without depth filter sheets.
- **Complete drainage** of filter due to unique filter plate design.
- **Safe operation:** No drip loss due to sealed external chamber and automatic hydraulic pressure regulation.
- **Easy handling** and **good cleanability** thanks to ergonomic design of chassis, and open support grid of filter plates; cake discharge downwards into tray underneath of filter package.
- **High flexibility:** The filter can be used as a sheet filter or as a pre-coat filter. For pre-coating/removing of high particle load, cake frames in different widths are available.
- BECO INTEGRA PLATE 600 EC filter is designed and manufactured according to the **cGMP guidelines** and the **FDA regulations**.

Customization or additional components allow for system expansion or complete systems according to customer specifications.



#### Technical Data

<b>Operating pressure:</b>	max. 87 psi (600 kPa/6 bar)
<b>Differential pressure:</b>	max. 58 psi (400 kPa/4 bar)
<b>Operating temperature:</b>	max. 284 °F (140 °C)
<b>Filter area per sheet:</b>	3.6 ft <sup>2</sup> (0.33 m <sup>2</sup> )

#### Materials

<b>Product contacting parts:</b>	Stainless steel AISI 316L/AISI 316Ti
<b>Other parts:</b>	Stainless steel AISI 304
<b>Gaskets/O-rings:</b>	Silicone, EPDM, Viton

Type	Filter area		Cake volume*		Chassis size	Filter plates/ cake frames	Depth filter sheets 23.0 x 26.8 in (584 x 680.5 mm)	Weight	
	ft <sup>2</sup>	(m <sup>2</sup> )	gal	(l)				qty	qty
<b>600 SF</b>	215.3	(20.00)	-	-	1	29/30	60	3,477	(1,577)
	387.5	(36.00)	-	-	2	53/54	108	4,306	(1,953)
	538.2	(50.00)	-	-	3	74/75	150	4,996	(2,266)
<b>600 ASF</b> with 0.31" (8 mm) cake frames	215.3	(20.00)	7.9	(30)	1	29/30	60	3,448	(1,564)
	387.5	(36.00)	14.3	(54)	2	53/54	108	4,259	(1,932)
	538.2	(50.00)	19.8	(75)	3	74/75	150	4,930	(2,236)
<b>600 ASF</b> with 0.71" (18 mm) cake frames	150.7	(14.00)	24.0	(91)	1	20/21	42	3,404	(1,544)
	272.7	(25.33)	43.0	(164)	2	37/38	76	4,182	(1,897)
	380.3	(35.33)	60.5	(229)	3	52/53	106	4,833	(2,192)
<b>600 ASF</b> with 0.98" (25 mm) cake frames	129.2	(12.00)	31.7	(120)	1	17/18	36	3,455	(1,567)
	229.6	(21.33)	56.3	(213)	2	31/32	64	4,246	(1,926)
	315.7	(29.33)	77.4	(293)	3	43/44	88	4,892	(2,219)
<b>600 ASF</b> with 1.57" (40 mm) cake frames	93.3	(8.67)	39.9	(151)	1	12/13	26	3,475	(1,576)
	165.0	(15.33)	70.8	(268)	2	22/23	46	4,268	(1,936)
	229.6	(21.33)	98.5	(373)	3	31/32	64	4,956	(2,248)

\* usable cake volume

#### Filter Plates/Cake Frames

	Width		Cake volume*		Weight	
	in	(mm)	gal	(l)	lb	(kg)
Filter plate	0.31	(8)	-	-	14.3	(6.5)
Feeding plate	0.31	(8)	-	-	14.3	(6.5)
Cake frame	0.31	(8)	0.26	(1.00)	13.5	(6.1)
Cake frame	0.71	(18)	1.14	(4.33)	23.3	(10.5)
Cake frame	0.98	(25)	1.76	(6.66)	32.2	(14.6)
Cake frame	1.57	(40)	3.08	(11.66)	51.6	(23.4)


\* usable cake volume

#### Connections

<b>Filter inlet:</b>	DN 65, DIN 11851/ISO 2 ½"
<b>Filter outlet:</b>	DN 65, DIN 11851/ISO 2 ½"
<b>External chamber:</b>	DN 15, DIN 11851/ISO ¾"

#### Other Designs

Flange according to DIN 2633/ISO or tri-clamp on request

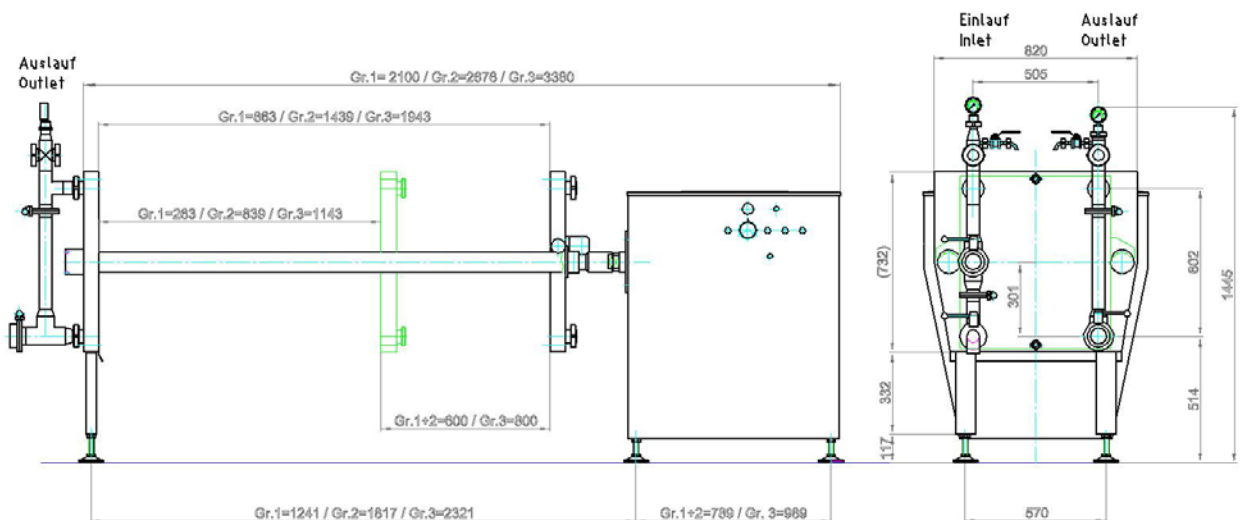



**North America**  
44 Apple Street  
Tinton Falls, NJ 07724  
Toll Free: 800 656-3344  
(North America only)  
Tel: +1 732 212-4700

**China**  
No. 3, Lane 280,  
Linhong Road  
Changning District, 200335  
Shanghai, P.R. China  
Tel: +86 21 5200-0099

**Europe/Africa/Middle East**  
Auf der Heide 2  
53947 Nettersheim, Germany  
Tel: +49 2486 809-0  
  
Friedensstraße 41  
68804 Altlußheim, Germany  
Tel: +49 6205 2094-0

**Singapore**  
4 Loyang Lane #04-01/02  
Singapore 508914  
Tel: +65 6825-1668

An den Nahewiesen 24  
55450 Langenlonsheim, Germany  
Tel: +49 6704 204-0

**Brazil**  
Av. Julia Gaioli, 474 –  
Bonsucesso  
07251-500 – Guarulhos, Brazil  
Tel: +55 11 2465-8822

**For more information, please  
email us at [filtration@eaton.com](mailto:filtration@eaton.com)  
or visit [www.eaton.com/filtration](http://www.eaton.com/filtration)**

EN  
1 A 2.6.11.2  
08-2015

© 2015 Eaton. All rights reserved. All trademarks and registered trademarks are the property of their respective owners. All information and recommendations appearing in this brochure concerning the use of products described herein are based on tests believed to be reliable. However, it is the user's responsibility to determine the suitability for his own use of such products. Since the actual use by others is beyond our control, no guarantee, expressed or implied, is made by Eaton as to the effects of such use or the results to be obtained. Eaton assumes no liability arising out of the use by others of such products. Nor is the information herein to be construed as absolutely complete, since additional information may be necessary or desirable when particular or exceptional conditions or circumstances exist or because of applicable laws or government regulations.